


Triple Rinse Procedure

- 1** Fill bottle with water to 10% of its capacity.

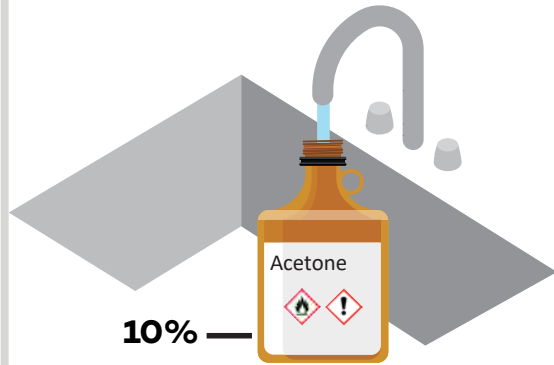


- 2** Replace cap, then shake and swirl the bottle.



- 3** Pour rinse water into a labeled waste container.



- 4** Repeat rinse process two more times.

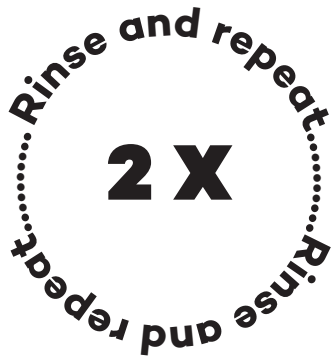

- 5** Deface original label and add Triple Rinse label.



- 6** Place bottle in hallway for custodians to recycle.



Do not triple rinse containers with sludge, solidified material, more than residual chemicals or septum caps. Submit as chemical waste instead.

Scan the QR code to watch the Triple Rinse video.

